





LABOMED, INC.

www.labomed.com
spectro@labomed.com

LB-1250 LCD Digital Biological Microscope Accessories

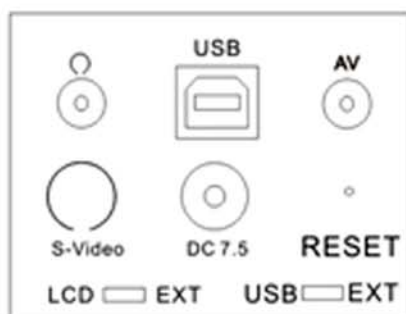

	Audio Jack		USB Output		RCA Video Output
	S-Video Output		Power Supply Input		Reset System
LCD EXT	LCD/Video Output	USB EXT	USB Output Switch		

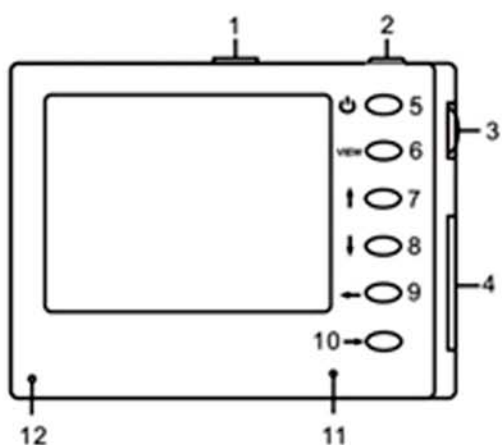

1. Screen latch release button, releases the screen when it is closed
2. Capture button, takes a still image, starts or stops capturing video
3. Menu/zoom direction button, press for menu/ok, push up/down for zoom in/out
4. SD card slot, external memory expansion slot
5. Power button, turns on and off the camera and LCD screen
6. View button, switches between live video and stored image viewer
- 7、8、9、10 Direction button, control the menu or image in viewer mode
11. Indicator light, blink red to acknowledge menu changes
12. MIC (Microphone), play the audio recorded